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OPERATING INSTRUCTIONS

1. SUPPLIED WITH	7
2. SECURITY ADVICE.....	7
3. TAKING INTO OPERATION.....	8
4. FUNCTIONALITY	8
5. SERVICE / MAINTENANCE	8
6. DISTURBANCES / REPAIR	8
7. WASTE DISPOSAL	9
8. SPECIFICATIONS	9
9. CUSTOMIZATIONS.....	9

SICCO Extractor Hood
Cat.No.: V 1960-02

SICCO Extractor Hood
Cat.No.: V 1960-04

Please read the information listed below thoroughly. We urgently ask you to observe the instructions about security, use and maintenance.

The SICCO Extraction Hood allows production of chemical compounds and provides safe protection from contaminating substances. The Extraction Hood can easily be integrated into your existing exhaust air system on site. All panels are made of high quality, transparent polycarbonate panels, except the top panel is made of polypropylene. The aluminium profiles are joint together with connectors made of polyamide.

1. Supplied with

Extractor Hood	Operating Instructions
1	1

2. Security advice

- » Do not take a damaged Extractor Hood into operation.
- » Provide a safe stand for the Extractor Hood.
- » SICCO Extractor Hood shall be taken into operation only in rooms with sufficient ventilation.
- » The ambient temperature range of $-20\text{ }^{\circ}\text{C}$ to $+70\text{ }^{\circ}\text{C}$ shall not be exceeded.
- » Do not expose the Extractor Hood to direct sunlight.
- » Place the Extractor Hood in a reachable working height.
- » Check the correct connection to your exhaust air system.
- » Check whether the performance of your exhaust air system is sufficient for the use of the Extractor Hood.
- » Do not lean the upper part of your body into the interior of the Extractor Hood.
- » Do not put the aspirator hose under tensile or compressive stress.
- » Please check after any absences whether the security instructions are still observed.
- » Do not work with any gas which is heavier than the ambient air.
- » This operating instruction cannot reflect all possible dangers which might occur during use. Users should check and observe all safety regulations.

SICCO **Extractor Hood**

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» Remove the packing and the protection cover and place the Extractor Hood on a horizontal, plain bench.

» Connect the Extractor Hood to your exhaust air system with suitable tube or tubing via the port on the top side. Secure the connection to avoid unintentional disconnection.

4. Functionality

» The Extractor Hood is connected to your existing exhaust air system by the connector port on its top side. Persons and the ambient atmosphere is protected from contamination by the transparent side walls.

5. Service / Maintenance

» Use only soft clothes for cleaning. Do not use any detergents which may contain solvents or which scrub.

6. Disturbances / Repair

» The following information shall help you to recognise and, if possible, to eliminate possible reasons for disturbances by yourself. If you are not able to handle the disturbance, please give us a call. We will then try to help you this way. If then, a solution to your problem cannot be found, we will ask you to return the cleaned glove box in its original packing to us. Then we will check, whether a cost-effective repair is possible.

Objection	Reason and help
The aspiration performance is too low.	» Foreign bodies can be stuck in the connector port or the aspirator tube. Remove them. » The distance from the chosen work place to the exhaust air system is too big. Choose an alternative place which is nearer to the exhaust air system. » The chosen work place can be situated at a minor line of your exhaust air system. Choose an alternative place and connect the extractor hood to the major line.

7. Waste Disposal

The Extractor Hood as well as its packing are made of recyclable materials. Please take further information on the materials used from the following specifications:

8. Specifications

	Extractor Hood Cat.No.: V 1960-02	Extractor Hood Cat.No.: V 1960-04
Overall dimensions: W x H x D	900 x 965 x 600 mm	1200 x 965 x 600 mm
Usable interior space: W x H x D	860 x 880 x 560 mm	1160 x 880 x 560 mm
Opening: W x H	545 x 600 mm	850 x 600 mm
Weight	20 kg	23 kg
Connector Ø	DN 160	
Material frame	aluminium with connectors made of polyamide	
Material panels	PC (Polycarbonate)	
Material top	PP (Polypropylene)	

9. Customizations

If you need an Extractor Hood adapted to your special needs, just send us an enquiry stating your requirements and dimensions. As manufacturer we are able to offer customized versions.

In general, the following modifications of our standard Extractor Hoods are possible:

- » Adaptation of the connector on the top side (without additional cost).
- » Additional connectors for ventilation or aspiration
- » Custom outer dimensions.
- » Panels in alternative materials
- » Antistatic version.

Please contact us for any questions regarding a customization.

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BOHLENDER GmbH