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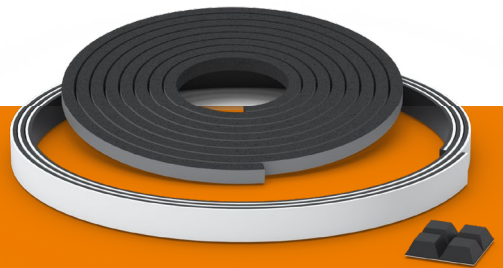
225 g

## OPERATING INSTRUCTIONS

1. SUPPLIED WITH .....	12
2. SECURITY ADVICE .....	12
3. FUNCTIONALITY .....	12
4. TAKING INTO OPERATION .....	13
5. SERVICE / MAINTENANCE .....	18
6. DISTURBANCES / REPAIRS .....	18
7. WASTE DISPOSAL .....	19
8. SPECIFICATIONS .....	19

**SICCO Maintenance Kit  
(Star-Desiccators)**  
Cat.No.: V 1778-02

**SICCO Maintenance Kit  
(Big Star and Maxi-Desiccators)**  
Cat.No.: V 1779-02

**Please read the information listed below thoroughly. We urgently ask you to observe the instructions about security, use and maintenance.**

By using the SICCO Gas-saving valve you can reduce your gas consumption significantly. The aeration tubing is closed nearly gas-tight. During gas-filling the gas-saving valve automatically opens due to the gas flow to prevent the accrument of overpressure inside your glove box.

## 1. Supplied with

### Maintenance Kit for Star-Desiccators

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1 Operating instructions

## 2. Security advice

- » Provide a safe stand for the desiccator during maintenance operation.
- » Run sharp tools always away from the body.
- » Use detergents in well-ventilated rooms only.

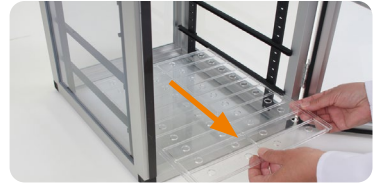
## 3. Functionality

The door is securely closed by means of magnetic tapes. The inside of the desiccator is reliably sealed against influences from the outside by means of sealing tape. The self-adhesive rubber feet provide a safe stand for your desiccator.


## SICCO Maintenance Kit

### 4. Taking into operation

- » Remove all components and the silica gel from the desiccator.
- » Remove all shelves from the desiccator.

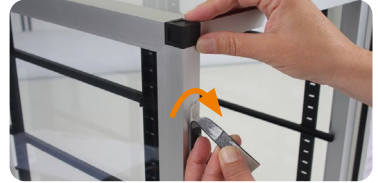- » Pull off the magnetic tape from the door and from the housing of the desiccator.

- » Only in rare cases the magnetic tape can be pulled off residue-free. There are two possible ways to remove the adhesive tape.
- » Peel off the adhesive tape with your thumb lengthwise.

- » Or use a knife to cut off the adhesive tape from the frame.
- » **Attention:** Cut away from your body. To get a good result the angle of the knife during cutting is crucial.

- » Remove the sealing tape accordingly.

- » Remove all adhesive residues. Use only soft cloths for cleaning. Do not use any detergents which may contain solvents or which scrub. You can use all common neutral detergents (pH 7).

## SICCO Maintenance Kit

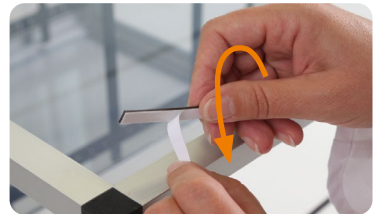
- » Mark the length of the aluminium profiles on the magnetic tapes.
- » Depending on the version the lengths are as follows:
  - Star- and Star Horizontal: one piece of type A and one piece of type B length 480 mm each.
  - Tower-Star: one piece of type A and one piece of type B length 690 mm each.
  - Big-Star: Two pieces with a length of 500 mm each.
  - Maxi 2: Four pieces with a length of 500 mm each.
  - Maxi 1: Two pieces with a length of 1030 mm each.


- » Cut the magnetic tapes to the required length by using scissors.

- » Pull off the protection film of the magnetic tape for approx. 10 cm.

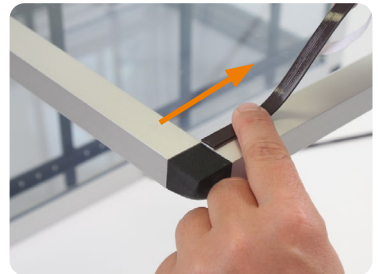- » Stick the magnetic tape flush with the inner edge of the aluminium profile of the housing. Pull the protection film further off. For versions Star, Star-Horizontal and Tower-Star the direction of the magnetic tape is random.

- » **Attention:** For versions Big-Star-, Maxi 1- and Maxi 2 it is crucial to stick the magnetic tape in the right direction.
- » Mount one magnetic tape to the housing of the desiccator. Take now the magnetic tape which you want to stick to the door and lay it on top of the already assembled magnetic tape.

## SICCO Maintenance Kit

- » If the direction of the second magnetic tape is wrong it will be offset.

- » Turn the magnetic tape top down.
- » If the direction is correct it will lay flush with the already mounted magnetic tape.

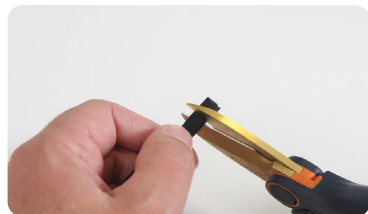
- » Mark the upper position with a suitable pen.

- » For versions Big-Star- and Maxi stick the marked magnetic tape on the door. Make sure that the marked part is at the top.
- » Stick the magnetic tape flush with the inner edge of the door on the aluminium profile. For versions Star-, Star-Horizontal- and Tower-Star the direction of the magnetic tape is not relevant.

- » If required cut one end of the sealing tape rectangular.

## SICCO Maintenance Kit

- » Remove approx. 10 cm of the protective film of the adhesive side of the sealing tape.
- » Start in the center of the hinge and stick the sealing tape on the frame. Keep a distance of approx. 1 mm to the outer edge of the aluminium profile.

- » Stick the sealing tape on the door. Be careful not to stretch or compress the sealing tape.

- » Stick the sealing tape round the corner. Make sure that the curved part of the sealing tape is placed on the corner connection.

- » Do not remove the last 10 cm of the protective film of the adhesive side of the sealing tape.
- » Overlap the beginning and the end of the sealing tape. Make sure to align the end of the sealing tape with the beginning before cutting.

- » Remove the protective film of the adhesive side and stick both ends of the sealing tape seamless together.

## SICCO Maintenance Kit

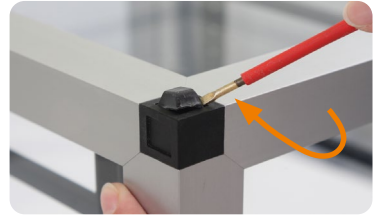
- » Close the door and visually check if the sealing tape is positioned correctly. Make sure to check all four sides.


- » **Tip:** Stick a small piece of paper between the sealing tape of the door and the frame of the desiccator. Pull the paper along the sealing tape. If you can feel a resistance on all four sides of the door, the desiccator is tight.


- » For versions of the Star-series remove the old rubber feet of the desiccator, e.g. by using a screwdriver with a 3,0 mm blade. Always run the screwdriver away from your body.

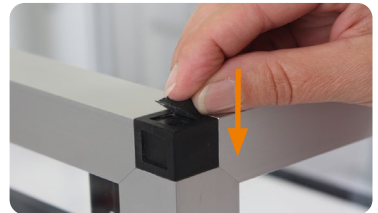- » Remove the protective film from the new rubber feet and stick them in the cavity of the corner connections. Press on the rubber feet.

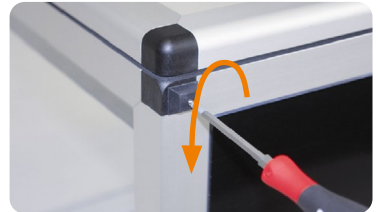- » For versions of the Big-Star-series remove all fastening screws of the rubber feet by turning them counterclockwise. Use a crosstip screwdriver size 1.
- » Keep the fastening screws for reuse.


## SICCO Maintenance Kit

- » Cut off the old rubber feet with a knife.

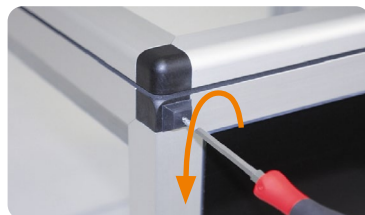
- » Stick the new rubber feet in center on the corner connections. Additionally use the fastening screws to fix the rubber feet. Screw them in by turning them clockwise. The screws may not project.

- » For versions of the Maxi-series remove the casters by drilling the blind rivet of the casters with a drill bit  $\varnothing$  5 mm until the rivet head gets loose. Attention: Risk of injury! If necessary ask an expert to execute this step.

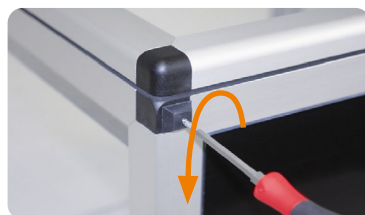- » Stick the rubber feet in center on the corner connections. Additionally use a countersunk with a  $\varnothing$  2,5 mm and a length of approx. 15 mm to fix each of the rubber feet. The screws are not included in delivery, a pre-drill is not necessary for this step.

### 5. Service / Maintenance

Use only soft clothes for cleaning. Do not use any detergents which may contain solvents or which scrub. You can use all common neutral detergents (pH 7).

### 6. Disturbances / Repairs

The following information shall help you to recognise and, if possible, to eliminate possible reasons for disturbances by yourself. If you are not able to handle the disturbance, please give us a call. We will then try to help you this way. If then, a solution to your problem cannot be found, we will ask you to return the cleaned desiccator in its original packing to us. Then we will check whether a cost-effective repair is possible.

# SICCO Maintenance Kit

## Objection

## Reason and help

The door cannot be closed.

» The magnetic tape was not stuck correctly. Check the position of both magnetic tapes according to the steps described under point 4. The magnetic tapes have to be stuck on the inner edge of the aluminium profiles.

**Info:** The magnetic tapes can be used only once.

» The sealing tape was not stuck correctly. Check the position of the sealing tape according to the steps described under point 4. The sealing tape has to be stuck on the outer edge of the door. **Info:** The sealing tape can be used only once.

The door cannot be closed tightly.

» The sealing tape has been stretched while sticking it on the door. A stretched sealing loses its elasticity and therefore cannot seal correctly. Exchange the sealing tape.

**Info:** The sealing tape can be used only once.

» There is another reason which cannot be handled with a Maintenance Kit for Desiccators. Check the door and the hinge-joints for deformation.

The desiccator jiggles.

» The rubber feet are not glued tightly. If you have a Star-Desiccator, check if all four rubber feet are glued into the slots. Remove the wrongly glued feet and glue them correctly into the slot. Use an appropriate glue such as Epple-Bond 8212. If you have a Big-Star-Desiccator, check if the fixing screws overlap. If they do, screw them in clockwise.

» There is another reason which cannot be handled with a Maintenance Kit for Desiccators. Check if your base is flat.

## 7. Waste Disposal

The Maintenance Kit for Desiccators as well as the packing consist of recyclable materials. Please take further information about the materials used from the following specifications.

## 8. Specifications

### SICCO Maintenance Kit for Desiccators

Cat.No.:	V 1778-02	V 1779-02
Use:	Star-Desiccators	Big-Star and Maxi Desiccators
Magnetic tape, self-adhesive:	0,75 m Type A, Width=8 mm 0,75 m Type A, Width=8 mm	2,5 m, Width=13 mm
Sealing Tape, self-adhesive:	2,0 m	5,0 m
Rubber feet, self-adhesive:	4 Pieces	4 Pieces
Material:	Cellular rubber, Polyurethane, Isoflex	
Weight:	125 g	225 g

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# Special Requirements? Custom-Made

## You are looking for something very particular? Something that even our giant portfolio of thought-through solutions does not cover?

No problem. As developer and producer we offer you the possibility of an individual production, as desired. This can be made faster and more economical than expected.

Just talk to our specialists about your ideas - we advise and support you already beginning with the construction and produce than material-appropriate according to your specifications. And this already from numer of pieces 1 at an unexpected price than you imagine.

For this, we only need a drawing (a rough sketch is sufficient) and some information.

## The check list for your tailor-made dessiccator/ glove box:

- ✔ What is the name of your dessiccator/ glove box?
- ✔ For which application do you need the desiccator/ glove box?
- ✔ Which outer dimensions do you need?
- ✔ What do you want to store?
- ✔ Which panel material do you require?
- ✔ What kind of shelves do you require?
- ✔ What will be the maximum weight load per shelf?
- ✔ What kind of accessories do you require (cylinder look, cable lead-ins, etc.)?
- ✔ Which number of pieces do you require?
- ✔ What is the budgetary price range per desiccator that should not be exceeded?

**Do you have a special request?**  
Just call us: **+49 (0) 93 46-9286-0**

Or just send us a drawing (rough scetch is sufficient) and some information to **info@sicco.de**. We will come back to you to discuss the details and establish a free offer.

Also made by Bohlander.

Best performance for your lab.

# Professional High-Performance Fluoroplastic Labware

PTFE | PFA | FEP

Stirrer Shafts, Magnetic StIRRung Bars or Tubing are only three of more than 1000 innovative BOLA products made of high performance plastics. For efficient, precise and safe work in your lab.

Ask for your catalogue or get more information:  
[www.bola.de/en](http://www.bola.de/en) or +49 (0) 93 46-92 86-0

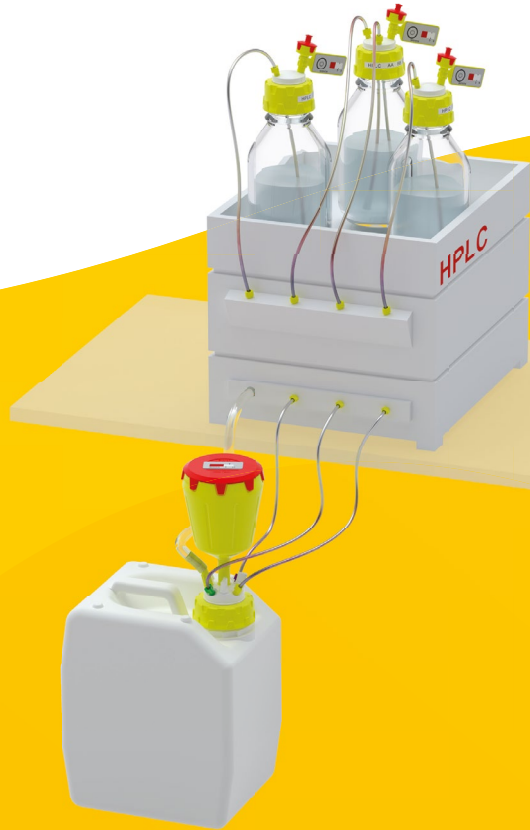
A PRODUCT BRAND OF BOHLENDER

[bola.de/en](http://bola.de/en)

Also made by Bohlender.

# b.safe

## Sophisticated Safety Solutions for HPLC.

**Easy. Better. Quicker. Directly from the manufacturer.**

When it comes to quality and safety, we leave nothing to chance: we develop, design and manufacture HPLC solutions in-house.

Ask for your catalogue or get more information:

**[www.bsafede/en](http://www.bsafede/en) or +49 (0) 93 46-92 86-0**

# SICCO

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**BOHLENDER GmbH**

Waltersberg 8

D 97947 Grünsfeld

+49 (0) 93 46-92 86-0

[info@sicco.de](mailto:info@sicco.de)

[www.sicco.de](http://www.sicco.de)