Bedienungsanleitung
Operating Instructions
Mode d'emploi

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
BOLA SAMPLING PUMP, Cat.-No.: A 124-16 A 124-16

Operating Instructions

1. Supplied with	5
2. Security Instructions	5
3. Taking into operation	5
4. Mode of function	5
5. Service / Maintenance	5
6. Disturbances / Repairs	5
7. Waste Disposal	6
8. Specifications	6
9. Spare Parts / Accessories	6

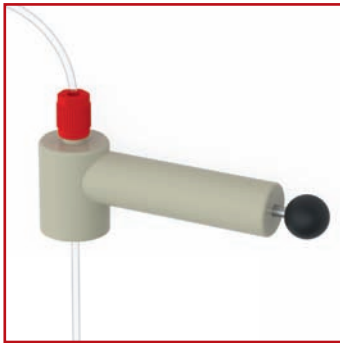Please read the information listed below thoroughly. We urgently ask you to observe the information about security, use and maintenance. BOLA Sampling Pumps are suitable for drawing and filling of liquids also at locations difficult to access. They can also be used for high viscous liquids such as oils.

BOLA Sampling Pump

1. Supplied with

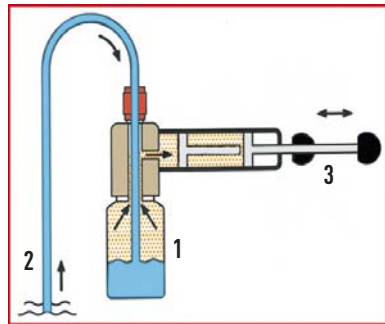
- 1 BOLA Sampling Pump
- 1 BOLA Laboratory Screw Joint for thread GL 18 and tubing O. D. 6 mm
- 2 Meters PTFE Tubing I. D. 4 x O. D. 6 mm
- 1 Operating Instructions

2. Security Instructions

- » Do not take a damaged pump into operation.
- » Wear protective gloves for pumping aggressive liquids.

3. Taking into operation

- » Screw the sampling pump onto your bottle (1) by turning it clockwise.
- » At least approx. 2 cm of the tubing have to extend in the bottle.
- » Immerse the tubing in the liquid (2) that has to be pumped.
- » For pumping, pull the handle (3) several times.
- » Approx. 50 ml per stroke are transported.

4. Mode of function

By pulling the handle (3), a slight vacuum is produced in the bottle. This vacuum pulls the sample (2) into the bottle (1). The sample is exposed only to the PTFE tubing and the collection vessel. Any glass, PTFE or PFA bottle with a thread GL 45 and a capacity from 100 ml up to 2000 ml can be used as collection vessel.

5. Service / Maintenance

Purge the pump and the tubing with a suitable cleaning liquid (e. g. clear water) after use.

4. Disturbances / Repairs

The following information shall help you to recognise and, if possible, to eliminate possible reasons for disturbances by yourself. If you are not able to handle the disturbance, please give us a call. We will then try to help you this way. If then, a solution to your problem cannot be found, we will ask you to return the cleaned sampling pump in the original packing to us. Then we will check whether a cost-effective repair is possible.

BOLA Sampling Pump

Objection

The pump is not pumping.

Reason and help

- » The tubing on the pumping side is not immersed in the productt. Put the tubing end into the liquid.
- » You have not made enough strokes. Pull again the handle for several times.
- » The mounted tubing is too long. Cut the tubing.
- » The tubing is blocked. Rinse the tubing with hot water.

7. Waste Disposal

The BOLA sampling pump as well as its packing consist of recyclable material.

Please take further information on the materials used from the following specifications:

8. Specifications

Bola Sampling Pump

Cat.-No.	Volume per stroke	Temperature resistance	Material	Bottle thread	Connecting thread for tubing	Tubing O. D.
A 124-16	ca. 50 ml	- 10 °C up to + 90 °C	Polyproylene stainless steel PTFE	GL 45	GL 18	6,0 mm

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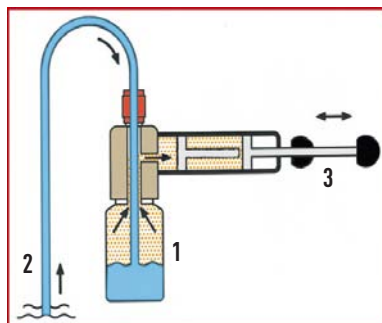
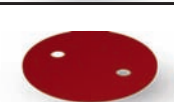
SPECIAL REQUIREMENTS? CUSTOMIZED!

You are looking for something very special?
Something that even our huge portfolio of
sophisticated lab solutions does not cover?

No problem:

As developer and producer, we offer the possibility
to produce individually according to your
requirement. This is faster, simpler and often
more economic than you can imagine. Just talk
to our experts about your ideas – we advise you
and support you already during the construction
and produce suitable for the material exactly
according to your specification. And this already
from quantity 1.

For this, we just need a drawing (a rough sketch is
sufficient) and some information.

Checklist for your customised product:

- » What is the article name?
- » In which application should the article be used?
- » What dimensions should the article have?
- » Are there any specific material specifications?
- » In which temperature range should the
article be used?
- » What chemical stresses is the article exposed to?
- » In which quantities is the article required?
- » What cost per piece should the article not exceed?

You have a special request?
Call us on: **+49 (0) 93 46-92 86-0**

Or just send us a drawing (a rough sketch is
sufficient) and some information to
info@bola.de. We will then contact you to
discuss further details and establish an offer
for you free of charge.

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BOHLENDER GmbH