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SICCO **Auto-Desiccator Wall**
Cat.No.: V 1995-08

SICCO **Auto Glove Box with Transfer Chamber** Cat.No.: V 1755-08

Please read the information listed below thoroughly. We urgently ask you to observe the instructions about security, use and maintenance.

SICCO Automatic Desiccators are made for long-term storage of reference samples, electronic components, photo equipment, documents , valuable assets and many more. The automatic drying unit reduces the relative humidity inside the desiccator to below 20 %. The drying units work maintenance-free. The panels of SICCO Auto-Desiccators are made of high-quality and transparent plastic material respectively of Borosilicat glass 3.3 for model Vitrum with a strong aluminium frame and connectors made of polyamide.

1. Supplied with	Shelves made of acrylic glass	Shelves made of stainless steel	Rails made of polyamide	Rails made of aluminium	Hygrometer	Operating instructions
Auto-Star-Desiccator	4		8		1	1
Auto-Star-Desiccator Vitrum		4		8	1	1
Auto-Big-Star-Desiccator		2	4		1	1
Auto-Maxi-1-Desiccator		4	8		1	1
Auto-Maxi-2-Desiccator		4	8		2	1
Auto-Desiccator-Wall		6	12		6	1
	Gloves size 9	The desiccant cartridges		Hygrometer	Operating instructions	
Auto Glove Box with Transfer Chamber	1x	3x		2x	2x	

2. Security Advice

- » Do not take a damaged desiccator into operation.
- » SICCO Automatic-Desiccators are suitable only for 100-230 V AC.
- » Do not open the control unit.
- » Do not store any solvents or acids in the desiccator.
- » Do not put any explosive or inflammable products into the desiccator.
- » SICCO Desiccators are not suitable for overpressure or vacuum.
- » Provide a safe stand for the desiccator.
- » Place the desiccator in a reachable working height.
- » The ambient temperature range of -20 °C to + 70 °C respectively of -20 °C to + 120 °C shall not be exceeded.
- » SICCO Automatic-Desiccators are to be used in closed and dry rooms only.
- » Protect the Automatic-Desiccators from splash water.
- » Do not expose the desiccator to direct sunlight.
- » Remove only one shelf at once.

3. Taking into operation

» Remove the protection cover.

» For versions with rails made of aluminium you need to remove the protection film from the rails.

» Remove the protection film from both sides of each shelf.

» Opening the door: Hold the desiccator housing with one hand while you open the door by pulling the door grip with the other hand.

» **Please note:** Due to the strong magnetic lock, it is harder to open the door the longer it was closed before. This is normal and can be compensated by pulling stronger than in the everyday use.

» Hook the rails made of aluminium (2) into the lateral mounted suspension arrangements (snap-in pin (1) up wards)

» Mount the rails (2) horizontally in pairs. Pay attention to mount them in the same height. A slight rotation of the rail helps to fix the snap-in pin safely in the suspension arrangement (3).

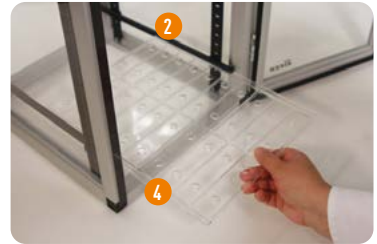
SICCO Auto-Desiccators

- » For models with rail (2) and suspension arrangement (3) made of plastic: Mount the rails (2) horizontally in pairs on both sides of the suspension arrangement (3). The engagement of the rail (2) is clearly audible. For a better orientation, the levels are marked by numbers on the suspension arrangement (3).

- » The shelves (4) can then be slid in on the rails (2). Mount another pair of rails (2) and slide in another shelf (4).
Please note: Mount the rails in pairs and slide in immediately a shelf to avoid that other rails can be fall down accidentally.

- » Open the battery case on the back of the hygrometer by pushing the cover downwards. Remove the battery safety strip and reinsert the battery. Close the battery case.

- » For desiccators with only one drying unit, connect the small plug (4) to the socket on the drying unit (6). Plug the power supply (5) into a 100-230 V socket.

- » For desiccators with several drying units, only the Schuko plug (8) needs to be plugged into a 100-230 V socket.

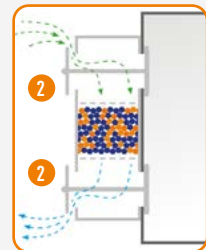
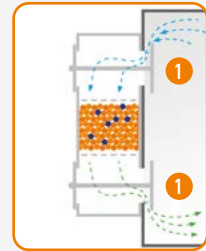
As soon as the drying unit is supplied with power, the valve slides open and close once, the fan starts to work and the status LED lights up green.

4. Functionality

The drying process is controlled by the automatic drying unit. It regulates the switch between drying the interior and the regeneration of the silica gel inside the drying unit. As soon as the drying unit is supplied with power, the green LED lights up and the fan runs continuously. Both valves open and close once, then stay in their default position where they are opened to the inside. (of the desiccator). The automatic cycle for dehumidification of the air inside the desiccator and regeneration of the desiccant takes about 40 minutes and operates continuously as follows:

- » During the 30 minute long drying cycle, both valves are in their default position (1). This means that the valves, and thus the drying unit, are open to the interior of the desiccator and closed to the outside. The fan draws air from the interior of the desiccator, feeds it through the desiccant and returns the dehumidified air to the desiccator.
- » Now the regeneration cycle for the regeneration of the desiccant starts. The green LED blinks during the regeneration cycle. The valves change their position (2) : they close the drying unit to the interior of the desiccator and open it towards the outside. The fan draws ambient air for 10 minutes and feeds it through the desiccant back again to the outside. Within the first eight minutes, the desiccant is heated up so that the absorbed humidity is released. The last two minutes of the regeneration cycle are used to cool down the desiccant to ambient temperature. Then the complete cycle starts from the beginning.

Functions of the digital Thermometer/Hygrometer:

Measuring range:	» Temperature – 10 °C to + 60 °C » Air humidity 20 % to 90 % relative humidity
Max. error in measurement:	» +/- 1 °C at –10 °C to + 50 °C » +/- 5 % relative humidity at 25 % to 75 % relative humidity and 0 °C to 50 °C » Attention: The mounted thermometer/hygrometer cannot display any values or data if the relative inside humidity of the desiccator is less than 25 % or higher than 75 %. To display exact values, we recommend using a thermometer/hygrometer of higher quality.
Battery:	» Please use a button cell LR44 or AG13.
Functionality of the Keys :	Indication of maximum / minimum values: » By pressing the MAX/MIN-key, the maximum value since the last reset will be displayed. » By pressing the MAX/MIN-key again, the minimum value since the last reset will be displayed. » The current values will be displayed after pressing MAX/MIN a third time. Reset of the MAX/MIN-memory: » Press CLEAR and the MAX/MIN-memory will be deleted. The current values will be stored instead.

Battery sign in the display:

- » If the battery is low, a battery flag is shown on the left bottom of the display. Exchange the battery to ensure best possible operation of the device.

5. Service / Maintenance

- » Use only soft clothes for cleaning. Do not use any detergents which may contain solvents or which scrub.
- » Keep the magnetic door strip clean.

6. Disturbances / Repairs

The following information shall help you to recognise and, if possible, to eliminate possible reasons for disturbances by yourself. If you are not able to handle the disturbance, please give us a call. We will then try to help you this way. If then a solution to your problem cannot be found, we will ask you to return the cleaned desiccator into the original packing to us. Then we will check, whether a cost-effective repair is possible.

Objection	Reason and help
Humidity inside the desiccator does not fall.	» The desiccator is not connected to the mains. Connect the plug to a socket (100-230 V AC). » The door is not closed tightly. Close the door. » The battery in the hygrometer is empty. Exchange the battery. » The built-in hygrometer might be faulty. Double-check the displayed values with an external hygrometer. Exchange the hygrometer.

7. Waste Disposal

The desiccators as well as their packing are made of recyclable materials. Please take further information on the materials used from the following specifications:

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SICCO Auto-Star-Desiccators

Cat.No.:

Drying Performance:

Power Consumption:

Mains Voltage:

Material panels:

Temperature resistance:

Material frame:

Lock:

Overall dimensions (W x H x D):

Excess length of drying unit:

Usable interior space (W x H x D):

Weight:

Max. number of shelves:

Max. all-over load per shelf:

Total all-over load:

V 1877-07

approx. 65 g moisture in 24 hours

max. 53 Watts

230 Volt AC

PMMA (Polymethylmethacrylate), rear panel: PC (Polycarbonate)

- 20° to + 70°C

Aluminium with connectors made of polyamide

Magnetic tape

310 x 525 x 375 mm

63 mm (back side)

260 x 480 x 330 mm

approx. 7,3 kg

26 pieces

10 kg

30 kg

SICCO Auto-Star-Desiccators Vitrum

Cat.No.:

Drying Performance:

Power Consumption:

Mains Voltage:

Material panels:

Temperature resistance:

Material frame:

Lock:

Overall dimensions (W x H x D):

Excess length of drying unit:

Usable interior space (W x H x D):

Weight:

Max. number of shelves:

Max. all-over load per shelf:

Total all-over load:

V 1823-07

approx. 65 g moisture in 24 hours

max. 53 Watts

230 Volt AC

Bororsilicate glass 3.3, rear panel: Aluminium

- 20° to + 120°C

Aluminium with connectors made of polyamide

Magnetic tape

310 x 525 x 375 mm

63 mm (back side)

260 x 480 x 330 mm

approx. 13,5 kg

26 pieces

10 kg

30 kg

SICCO Auto-Big-Star-Desiccators

Cat.No.:

Drying Performance:

Power Consumption:

Mains Voltage:

Material panels:

Temperature resistance:

Material frame:

Lock:

Overall dimensions (W x H x D):

Excess length of drying unit:

Usable interior space (W x H x D):

Weight:

Max. number of shelves:

Max. all-over load per shelf:

Total all-over load:

V 1894-07

approx. 106 g moisture in 24 hours per chamber

max. 40 Watts

230 Volt AC

PMMA (Polymethylmethacrylate), rear panel: PC (Polycarbonate)

- 20° to + 70°C

Aluminium with connectors made of polyamide

Magnetic tape

560 x 560 x 580 mm

63 mm (back side)

495 x 500 x 540 mm

approx. 22 kg

17 pieces per chamber

30 kg

80 kg

SICCO Auto-Maxi 1-Desiccators

Cat.No.:

Drying Performance:

Power Consumption:

Mains Voltage:

Material panels:

Temperature resistance:

Material frame:

Lock:

Overall dimensions (W x H x D):

Excess length of drying unit:

Usable interior space (W x H x D):

Weight:

Max. number of shelves:

Max. all-over load per shelf:

Total all-over load:

V 1914-01

approx. 220 g moisture in 24 hours

max. 212 Watts

230 Volt AC

PMMA (Polymethylmethacrylate), rear panel: PC (Polycarbonate)

- 20° to + 70°C

Aluminium with connectors made of polyamide

Magnetic tape

560 x 1150 x 580 mm

63 mm (back side)

495 x 1030 x 540 mm

approx. 31 kg

34 pieces

30 kg

160 kg

SICCO Auto-Maxi 2-Desiccators

Cat.No.:

Drying Performance:

Power Consumption:

Mains Voltage:

Material panels:

Temperature resistance:

Material frame:

Lock:

Overall dimensions (W x H x D):

Excess length of drying unit:

Usable interior space (W x H x D):

Weight:

Max. number of shelves:

Max. all-over load per shelf:

Total all-over load:

V 1916-01

approx. 115g moisture in 24 hours per chamber

max. 212 Watts

230 Volt AC

PMMA (Polymethylmethacrylate), rear panel: PC (Polycarbonate)

- 20° to + 70°C

Aluminium with connectors made of polyamide

Magnetic tape

560 x 1150 x 580 mm

63 mm (back side)

495 x 1030 x 540 mm per chamber

approx. 35 kg

17 pieces per chamber

30 kg per chamber

80 kg per chamber

SICCO Auto-Desiccators-Wall

Cat.No.:

Drying Performance:

Power Consumption:

Mains Voltage:

Material panels:

Temperature resistance:

Material frame:

Lock:

Overall dimensions (W x H x D):

Excess length of drying unit:

Usable interior space (W x H x D):

Weight:

Max. number of shelves:

Max. all-over load per shelf:

Total all-over load:

V 1995-08

approx. 115 g moisture in 24 hours per chamber

max. 636 Watts

230 Volt AC

PMMA (Polymethylmethacrylate), rear panel: PC (Polycarbonate)

- 20° to + 70°C

Aluminium with connectors made of polyamide

Magnetic tape

1080 x 1870 x 580 mm

100 (back side)

495 x 500 x 540 per chamber

approx. 100 kg

17 pieces per chamber

30 / 80 per chamber

80 kg per chamber

SICCO **Auto-Desiccators**

SICCO Auto Glove Box with Transfer Chamber

Cat.No.:

Drying Performance:

Power Consumption:

Mains Voltage:

Material panels:

Temperature resistance:

Material frame:

Lock:

Overall dimensions (W x H x D):

Excess length of drying unit:

Usable interior space (W x H x D):

Weight:

Total all-over load:

V 1755-08

approx. 160 g moisture in 24 hours

max. 212 Watts

230 Volt AC

PMMA (Polymethylmethacrylate), rear panel: PC (Polycarbonate)

- 20° to + 70°C

Aluminium with connectors made of polyamide

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Hereby declares that the aforementioned products comply with the requirements of the following directives:


Low Voltage Directive 2014/35/EU

The following directives were used for assessment:

ISO 12100: 2010
DIN EN ISO 12100
EN 61326-1 1997

The product fulfills the requirements provided that its intended use is ensured during installation and operation.

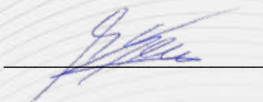Grünsfeld, 03.08.2026

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Desiccators and Drying Cabinets

Drying Cabinets, Desiccators and Glove Boxes

Stable | Functional | Custom-Made

For work in an isolated atmosphere and for dust- and moisture-free storage or storage with inert gas.

Ask for your catalogue or get more information:

www.sicco.de/en or +49 (0) 93 46-92 86-0

A PRODUCT BRAND OF BOHLENDER

sicco.de/en

Special Requirements? Custom-Made

You are looking for something very particular? Somthing that even our giant portfolio of thought-through solutions does not cover?

No problem. As developer and producer we offer you the pdsibility of an individual production, as desired. This can be made faster and more economical than expected.

Just talk to our specialists about your ideas - we advise and support you already beginning with the construction and produce than material-appropriate according to your specifications. And this already from numer of pieces 1 at an inexpected price than you imagine.

For this, we only need a drawing (a rough sketch is sufficient) and some information.

The check list for your tailor-made dessiccator/ glove box:

- ✓ What is the name of your dessiccator/ glove box?
- ✓ For which application do you need the desiccator/ glove box?
- ✓ Which outer dimensions do you need?
- ✓ What do you want to store?
- ✓ Which panel material do you require?
- ✓ What kind of shelves do you require?
- ✓ What will be the maximum weight load per shelf?
- ✓ What kind of accessories do you require (cylinder look, cable lead-ins, etc.)?
- ✓ Which number of pieces do you require?
- ✓ What is the budgetary price range per desiccator that should not be exceeded?

Do you have a special request?
Just call us: **+49 (0) 93 46 92 86-0**

Or just send us a drawing (rough scetch is sufficient) and some information to **info@sicco.de**. We will come back to you to discuss the details and establish a free offer.

SICCO

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BOHLENDER GmbH

Waltersberg 8

D 97947 Grünsfeld

+49 (0) 93 46-92 86-0

info@sicco.de

www.sicco.de