**Bedienungsanleitung**  
**Operating Instructions**  
**Mode d'emploi**

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» siehe Seite 16

## **BOLA Magnetic Stirrer Heads**

**BOLA Magnetic Stirrer Heads with Ground Joint, Cat.-No.: C 450-16 and C 450-24.**

**BOLA Magnetic Stirrer Heads with Flange, Cat.-No.: C 454-24.**

### **OPERATING INSTRUCTIONS**

<b>1. Supplied with</b>	<b>7</b>
<b>2. Security instructions</b>	<b>7</b>
<b>3. Mode of function</b>	<b>7</b>
<b>4. Taking into operation</b>	<b>7</b>
<b>5. Service / Maintenance</b>	<b>10</b>
<b>6. Disturbances/ Repairs</b>	<b>10</b>
<b>7. Waste disposal</b>	<b>10</b>
<b>8. Specifications</b>	<b>10</b>
<b>10. Spare parts / Accessories</b>	<b>10</b>

BOLA Magnetic Stirrer Heads with Ground Joint,  
Cat. No.: C 450-16 and C 450-25..

BOLA Magnetic Stirrer Heads with Flange,  
Cat. No.: C 454-24.

Please read the information listed below thoroughly. We urgently ask you to observe the information about security and use.

BOLA Magnetic Stirrer Heads with Ground Joint and BOLA Magnetic Stirrer Heads with Flange are suitable for stirring under vacuum. They can either be fit in a ground joint size 29 or on a flange size 25. The stirrer heads can accept stirrer shafts with a diameter of 8 mm. Stirring is effected without abrasion.

The product which is stirred is only in contact with PTFE and PFA which have a very good chemical resistance.

# BOLA Magnetic Stirrer Heads


## 1. Supplied with

- 1 BOLA Magnetic Stirrer Head (1)
- 1 Assembling aid (2)
- 1 Adaptor for assembling in a chuck (3)
- 1 Operating instructions

## 2. Security instructions

- » Do not take a damaged magnetic stirrer head into operation.
- » Only use the magnetic stirrer head up to the maximum speed of 800 rpm.
- » Never use the magnetic stirrer head (B) without globe stirrer coupling (A), Cat. No. C 399-12, or a comparable coupling.
- » There must be a gap of approx. 2 mm (see picture on the right side) between magnetic stirrer head (B) and globe stirrer coupling (A) for compensating heat expansion.
- » Only wear tight clothes to avoid being drawn in by rotating components.
- » If you have long hair, please wear a hairnet.
- » In case of temperature fluctuation, the correct fit of the stirrer shaft must be checked carefully.
- » Please only use tight gloves.
- » Only use the magnetic stirrer head for products which are in a temperature range between - 200 °C to +250 °C.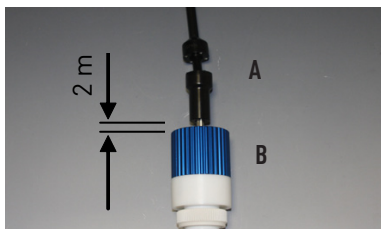

## 3. Mode of function

Due to the built-in permanent magnets in the blue drive shaft and in the rotor, the torque is transmitted from the motor to the stirrer shaft in the reactor. The casing of the blue drive shaft does not turn.

## 4. Taking into operation


- » Remove the nut (4) completely by turning it in counterclockwise direction. Hold the square drive journal with your other hand.


## BOLA Magnetic Stirrer Heads


» We recommend to use stirrer shafts with a clamping pin like shown on the picture.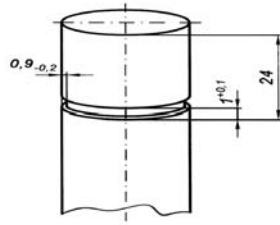


» If the stirring diameter of your stirrer shaft is bigger than the inner diameter of the ground joint or flange, the stirrer shaft must be introduced through the reactor lid from below.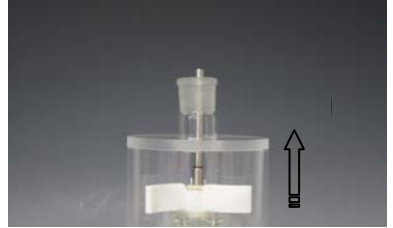


» Push the nut (4) over the stirrer shaft. The thread of the nut points to the end of the stirrer shaft.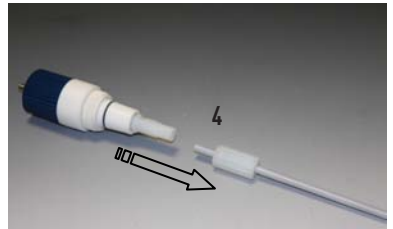

» First push the sealing ring (5) and then the clamp ring (6) on the stirrer shaft. The clamp ring must lock into the groove of the stirrer shaft. Observe the direction of assembly!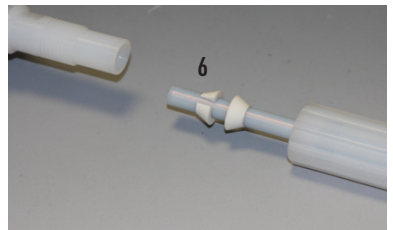

» If you have a short stirrer shaft, it is possible that the groove is inside the nut (4). In this case, take the assembling aid (2) and position the sealing ring (5) and clamp ring (6). The clamp ring must lock into the groove of the stirrer shaft.

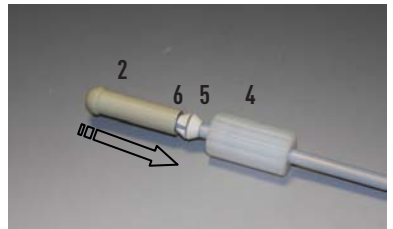» Push this component assembly carefully into the stirrer head. Tighten the nut (4) with two fingers on the shaft. Use a 14 mm flat spanner for locking.


## BOLA Magnetic Stirrer Heads

» Turn the pull nut (7) in clockwise direction completely to the top (only for ground joint version).


» Carefully put the stirrer head in the ground joint.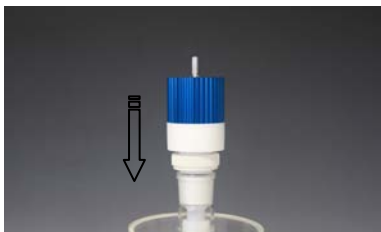

» For the flange version, use a suitable connector.

- » Assemble the stirrer coupling (A) between stirrer motor and stirrer head (B).
- » Adjust the speed control of the stirrer motor to 0 rpm. Approach the requested speed carefully by increasing slowly during stirring. Please observe the maximum speed of 800 rpm.

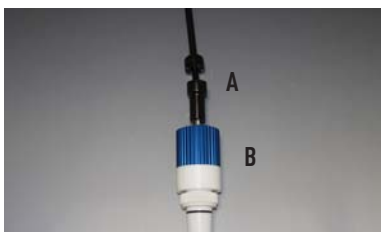» You can also fit the stirrer head in an 8 mm chuck. Please use the included adaptor (3) for this.

» A stuck stirrer head can be removed from the ground joint by turning the pull nut (7) in counterclockwise direction.

## BOLA Magnetic Stirrer Heads

### 5. Service / Maintenance

Please do not use abrasive detergents since these might damage the surface and result in a milkiness of the surface. You can use all common neutral detergents (pH 7). For a stronger contamination we recommend to use an alkaline detergent up to pH 12. The BOLA Magnetic Stirrer Head can be used without additional lubricants.

### 6. Disturbances/ Repairs

Due to their complex construction, we do not recommend you to disassemble the stirrer heads by your own, but to return it to us in cleaned condition. We will then check if a cost-effective repair is possible.

### 7. Waste disposal

The packing as well as the magnetic stirrer head are made of recyclable materials. Please take further details about the materials from the following specifications.

### 8. Specifications

#### BOLA Magnetic Stirrer Heads with Ground Joint

Cat. No.	Ground Joint	Dimensions	For stirrer shaft dia.	Material	Torque	Max. speed	Drive
C 450-16	NS 29/32	Ø 50 x 203 mm	8 mm	PPFA, PTFE Ceramic Aluminium Borosilicate glass Stainless steel 1.4305	20 Ncm	800 rpm	Outer square size 6  Adaptor Ø 8 mm
C 450-24		Ø 50 x 215 mm			40 Ncm		

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Cat. No.	Flange	Dimensions	For stirrer shaft dia.	Material	Torque	Max. speed	Drive
C 454-24	NW 25	Ø 92 x 215mm	8 mm	PPFA, PTFE Ceramic Aluminium Borosilicate glass Stainless steel 1.4305	60 Ncm	800rpm	Outer square size 6  Adaptor Ø 8 mm

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**CE Declaration of Conformity**

**EC-Declaration of Conformity in compliance with Machinery Directive 2006/42/EC**

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herewith declares that the following products:

**Magnetic Stirrer Heads (P-MRK), product group: C 520- ...**

**Magnetic Stirrer Heads (P-MRK) with Flat Flange, product group: C 522- ...**

**Magnetic Stirrer Heads with Ground Joint, product group: C 450- ...**

**Magnetic Stirrer Heads with Flange, product group: C 454- ...**

apply to the above mentioned directive including any amendments valid at the time this declaration was signed.

**The following directives have been applied:**

**Directive 2014/30 EU**

**Directive relating to electromagnetic compatibility,  
General requirements**

**The low voltage directive 2006/95/EC has been fulfilled according to 2006/42/EC terms of protection objectives.**

**The following harmonized standards have been applied:**

**DIN EN ISO 12100**

**Safety of machinery**

**ISO 12100:2010**

**Risk assessment and risk reduction**

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*As a manufacturing company we can make products to your specifications. Please send us a drawing or a sketch together with answers to the questions listed below.*

*Return this form with contact details by fax or e-mail and we will be in touch to discuss details and give you a free, no obligation quote.*

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BOHLENDER GmbH