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Operating Instructions

Mode d'emploi

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BOLA Fluoroplastic Spray, Cat.-No.: H 958-04

OPERATING INSTRUCTIONS

1. Supplied with	3
2. Hazard information	3
3. Safety information	4
4. Instruction for use	4
5. Mode of function	4
6. Waste disposal	4
7. Specifications	4

Please read the information listed below thoroughly. We urgently ask you to observe the information about security and use of this product.

1. Supplied with

- 1 Fluoroplastic Spray (400ml)
- 1 Operating Instructions

2. Hazard information

- » Extremely flammable aerosol.
- » Pressurized container: may explode if heated.
- » Causes serious eye irritation.
- » Harmful to aquatic life with long lasting effects.

BOLA Fluoroplastic-Spray

3. Safety information

- » Please do not take a damaged container into operation.
- » Keep out of the reach of children.
- » Keep away from heat/sparks/open flames/hot surfaces. No smoking.
- » Pressurized container: Do not pierce or burn, even after use.
- » Do not spray on an open flame or other ignition source.
- » Wear eye protection/face protection.
- » Do not breathe vapours/spray. Use only outdoors or in a well-ventilated area.
- » IF IN EYES: Rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing.
- » Protect from sunlight. Do not expose to temperatures exceeding 50 °C.
- » Do not pierce or burn even after use.
- » Keep away from any ignition source.

4. Instruction for use

- » Part must be clean, dry and free of grease, so that the film sticks on the surface.
- » Shake the can vigorously before every use.
- » Hold the can in direction of the part with a distance of approx. 20-25 cm. Spray several times if necessary.
- » If several layers shall be applied, each layer has to dry separately.
- » Polishing improves the anti-adhesive property.
- » The film is not scratch-proof. It can be removed by rubbing with a rough cloth or a brush

5. Mode of function

Fluoroplastic spray is used as dry, fat free separating agent and lubricant in industry, workshops and households. It is perfect for greasing parts, tools and devices (no freezing).

With fluoroplastic spray you obtain good sliding properties and a non-adhesive protective film. The sprayed-on film is heat resistant up to 260 °C. No smearing or soiling.

The fluoroplastic spray is CFC-free.

6. Waste disposal

- » The cans as well as their packing are made of recyclable materials. Dispose the content/container as per the local/regional/national/international regulations.

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BOHLENDER GmbH