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Operating Instructions

Mode d'emploi

» Mini-Stand-Hygrometer

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» Mini Hygròmetre

SICCO Mini-Stand-Hygrometer, Artikelnummer: V 1829-08

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SICCO Mini Hygrometer, Cat.-No.: V 1829-08

OPERATING INSTRUCTIONS

- | | |
|---|---|
| 1. Supplied with | 3 |
| 2. Taking into operation | 3 |
| 3. Functionality /Technical Information | 4 |

Please read the information listed below thoroughly. We urgently ask you to observe the instructions about use and maintenance.

1. Supplied with

- 1 SICCO Mini Hygrometer
- 2 Batteries 1,5V AAA
- 1 operating instructions

2. Taking into operation

» Take out the hygrometer and the batteries carefully.

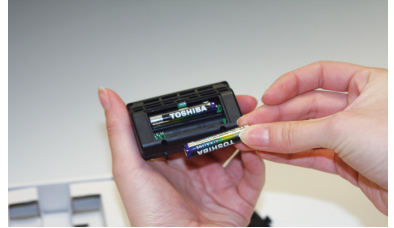
Sicco Mini Hygrometer


» Open the lid on the backside of the hygrometer.

» While inserting the batteries pay attention to the polarity. Corresponding symbols are inside the battery case.

» Click the lid back into place.

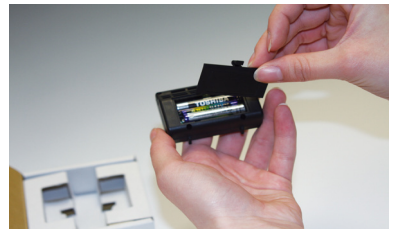3. Functionality /Technical Information

Measuring range:	» Temperature: 0 to 50 °C » Humidity 1 – 99 % rel. humidity
Max. measuring error:	» ± 1 °C at 25 °C » ± 3 % rel. humidity at 1 % to 80 % » The mounted hygrometer cannot display any values or data if the relative inside humidity of the desiccator is less than 1 % or higher than 80 %. To display exact values, we recommend using a hygrometer of higher quality.
Sensor drift:	» 0,2 °C in 4 years » 1 % rel. humidity in 4 years
Measuring rate	» 45 sec.
Resolution:	» 1 °C » 1 %
Battery indicator:	» The hygrometer has no battery indicator. With decreasing battery power the intensity of the LED-display is reduced. Therefore we recommend to check the hygrometer regularly and, if necessary, exchange the batteries.

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BOHLENDER GmbH