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Operating Instructions

Mode d'emploi

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OPERATING INSTRUCTIONS

1. SUPPLIED WITH	10
2. SECURITY INSTRUCTIONS	10
3. MODE OF FUNCTION	10
4. TAKING INTO OPERATION	11
5. SERVICE / MAINTENANCE	14
6. DISTURBANCES / REPAIRS	15
7. WASTE DISPOSAL	16
8. SPECIFICATIONS	16
9. SPARE PARTS / ACCESSORIES	16

Please read the information listed below thoroughly. We urgently ask you to observe the information about security and use.

BOLA Globe Stirrer Couplings are ideal for balancing misalignment of axes between agitator and stirrer shaft. They damp vibrations and increase the life expectancy of your stirrer bearing. BOLA Globe Stirrer Couplings are suitable for glass, metal or BOLA stirrer shafts.

BOLA Globe Stirrer Couplings

1. Supplied with

1 BOLA Globe Stirrer Coupling

For the following versions a reducing sleeve is included in delivery: cat.-no. C 394-02 / C 398-08 and cat.-no. C 394-03 / C 398-12

1 Operating instructions

2. Security instructions

- » Do not take a damaged Globe Stirrer Coupling into operation.
- » Make sure that the stirrer shaft is fixed tightly in the Globe Stirrer Coupling before starting a new stirring interval.
- » Make sure that the Globe Stirrer Coupling is fixed tightly in the agitator before starting a new stirring interval.
- » Only use the Globe Stirrer Couplings for the intended chucking diameters respectively for the intended form-fit supporting possibilities.
- » Do not use the Globe Stirrer Coupling without a stirrer bearing for safe guidance of the stirrer shaft.
- » Place the agitator in centre above the stirrer shaft.
- » No left-handed rotation is possible for the version with GL 10 thread. The stirrer shaft would be unscrewed if used with left-hand rotation.
- » Due to the indolence of the stirrer shaft and the turning medium you have to be careful that stirrer shafts with GL 10 thread are not unscrewed when you stop the agitator.
- » Start your stirring process by using the lowest possible rotation speed and slowly accelerate your agitator until the requested rotation speed is reached.
- » Slowly reduce the rotation speed to the lowest before you completely stop the agitator.
- » Only use the BOLA Globe Stirrer Couplings to the maximum permissible rotation speed of 1200 revolutions per minute.
- » Wear tight fitting clothing.
- » In case you have long hair we recommend to wear a hair net to avoid the danger of being pulled in.
- » If you wear gloves make sure they are tight fitting.
- » At least 30 mm of the pivot of the Globe Stirrer Coupling have to be clamped in the agitator.
- » Slide your stirrer shaft into the opening of the Globe Stirrer Coupling until it abuts on the ground.
- » For the version with a hexagonal reception shaft make sure to clamp on the hexagonal surfaces.
- » In case of temperature fluctuations make sure that the stirrer shaft fits tightly.
- » Never leave your equipment unsupervised during the stirring process.

These operating instructions cannot reflect all possible dangers which might occur during use. It's the users liability to observe security and health instructions and to fix corresponding restrictions before use.

3. Mode of function

A possible misalignment of axes between stirrer shaft and agitator can be balanced by means of a double ball-and-socket joint. Due to the material of the Globe Stirrer Couplings vibrations are damped. Depending on the version of the Globe Stirrer Coupling stirrer shafts can be accepted by either a collet chuck or a GL 10 thread. The version with inner square is ideal for a power transmission for example on a magnetic stirrer head.


BOLA Globe Stirrer Couplings

4. Taking into operation


- » Adjust the lowest possible rotation speed on the speed control.
- » Turn off the agitator.


- » For versions with a collet chuck open the clamping nut (1) by turning it counter-clockwise two times.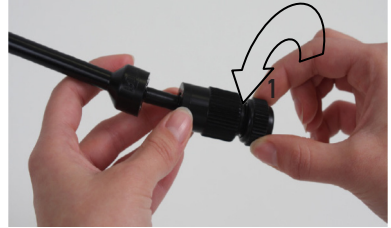

- » Please check if the chucking diameter of your stirrer shaft fits into the Globe Stirrer Coupling. Once the collet chuck is open the stirrer shaft should be easy to insert. Make sure that the shaft fits tightly without any gaps. Stirrer shafts with chucking diameter of 4 respectively of 10 mm can directly be inserted and fixed.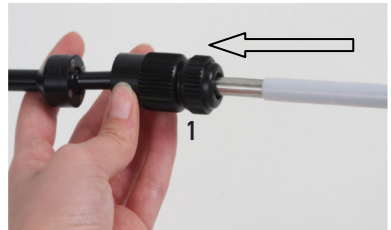

- » For stirrer shafts with a chucking diameter of 6,5 respectively of 8,0 mm a reducing sleeve (2) is included. If required slide the reducing sleeve (2) completely through the clamping (1) nut into the collet chuck.

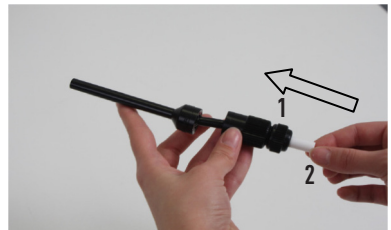- » The version with inner square, feed size 6mm, is ideal for applications with an appropriate stirrer shaft and it is designed to be used for example with a BOLA Magnetic Stirrer Head. The inner square of the Globe Stirrer Coupling should be easy to slide on the driving tap of your stirrer shaft or of your Magnetic Stirrer Head. The BOLA Magnetic Stirrer Head is available as an accessory, cat-no. C 520-... .




- » Due to the larger manufacturing tolerances of glass threads stirrer shafts with a GL 10 thread may be very easy to screw in. Fasten the stirrer shaft by turning it clockwise with two fingers.

- » **Attention:** The version with GL 10 thread is only suitable for right-handed rotation.



BOLA Globe Stirrer Couplings

- » These operating instructions cannot show all applications for the BOLA Globe Stirrer Coupling. The following assembly example shows the use with a stirrer shaft which has a chucking diameter of 8,0 mm. The stirrer shaft is guided by stirring bearing.
- » Station the agitator and your vessel on a stand.
- » Open the screw cap (3) of the stirrer bearing by turning it counter-clockwise two times. Slide your stirrer shaft through the stirrer bearing from below. The stirrer bearing is available as accessory, cat.-no. C 424-... .
- » Position the stirrer shaft inside the vessel on the requested level. Fasten the screw cap (3) by turning it clockwise with two fingers.
- » Generally the collet chuck of your agitator is hollow inside. Insert the Globe Stirrer Coupling completely into the collet chuck.
- » If necessary the pivot of the Globe Stirrer Coupling can be shortened by using a fine-toothed saw.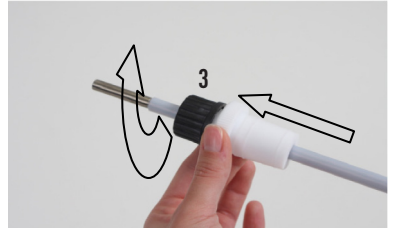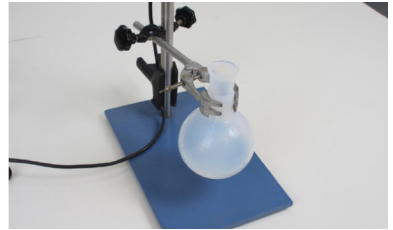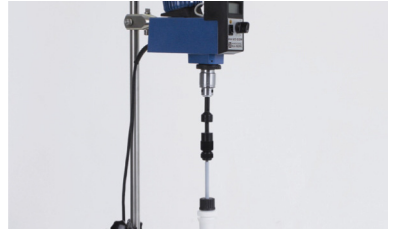

BOLA Globe Stirrer Couplings

- » Move your vessel with the mounted stirrer shaft and place it directly under the stirrer shaft.
- » Open the collet chuck of your agitator by turning it counter-clockwise two times.
- » Carefully pull the Globe Stirrer Coupling out of the collet chuck of your agitator and slide it on the stirrer shaft.
- » Fasten the collet chuck of your Globe Stirrer Coupling by turning the clamping nut (1) clockwise.
- » Close the collet chuck of the agitator by turning it clockwise. Use the belonging chuck key for this step.
- » Align the vessel with the mounted stirrer shaft and the Globe Stirrer Coupling as straight as possible.
- » Make sure that all fastening screws on the stand and on the agitator are fixed tightly.
- » Info: The BOLA Globe Stirrer Coupling can balance a maximum misalignment of axes of 10 mm.

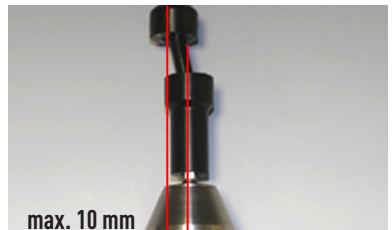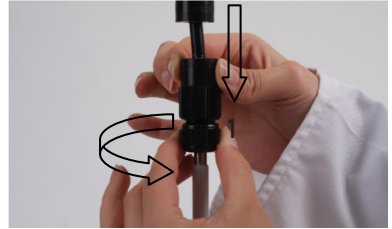BOLA Globe Stirrer Couplings

- » Adjust the lowest possible rotation speed on the speed control of the agitator and turn on the agitator. Slowly accelerate the agitator until the required rotation speed is reached.
- » **Attention:** The maximum speed is 1200 revolutions per minute.
- » Never leave the equipment unsupervised during the stirring process.
- » At the end of the stirring process slowly reduce the rotation speed to the lowest possible value. Then stop your agitator.

5. Service / Maintenance

Use only soft clothes for cleaning. You can use all common neutral cleaners (pH 7). In the case of a coarse contamination the clamping nut (1) and the reducing sleeve (3) can be removed and cleaned separately. The Globe Stirrer Coupling must not be immersed into the cleaning liquid completely. Otherwise the greasing inside which is maintenance free will be damaged.

6. Disturbances / Repairs

The following information shall help you to recognise and, if possible, to eliminate possible reasons for disturbances by yourself. If you are not able to handle the disturbance, please give us a call. We will then try to help you this way. If then, a solution to your problem cannot be found, we will ask you to return the cleaned Globe Stirrer Coupling to us. Then we will check whether a cost-effective repair is possible.

BOLA Globe Stirrer Couplings

Objection	Reason and Help
The stirrer shaft cannot be clamped in the collet chuck of the Globe Stirrer Coupling.	<ul style="list-style-type: none">» The chucking diameter of your stirrer shaft isn't compatible with the collet chuck. Use the included reducing sleeve. Depending on the version a reducing sleeve from Ø 10 to Ø 8 mm, respectively from Ø 10 to Ø 6,5 mm is included.» The reducing sleeve isn't compatible with your stirrer shaft. Use a stirrer shaft with a chucking diameter of 6,5 mm or 8,0 mm. Stirrer shafts with a chucking diameter of 4 and 10 mm are clamped without a reducing sleeve.» Residues inside the collet chuck prevent the correct clamping. Unscrew the clamping nut (1) by turning it counter-clockwise. Clean the collet chuck and the clamping nut (1) with a suitable detergent, for example with warm soap water.
The collet chuck cannot be opened.	<ul style="list-style-type: none">» Residues agglutinate the thread of the clamping nut (1). Immerge the Globe Stirrer Coupling with the clamping nut (1) ahead approx. 1 cm into hot soap water. Screw off the clamping nut (1) from the collet chuck by turning it counter-clockwise. We recommend to wear heat resistant gloves. <p>Attention: Risk of burns!</p>
The stirrer shaft does not turn.	<ul style="list-style-type: none">» The stirrer shaft is not clamped tight enough in the Globe Stirrer Coupling. Check whether the clamping nut (1) is sufficiently tight. If necessary fasten the clamping nut (1) by turning it clockwise with two fingers.» The stirrer bearing was closed too tight and thus prevents the stirrer shaft from turning. Loosen the screw cap of the stirrer bearing until the stirrer shaft can be turned.» The medium is too viscous. BOLA Globe Stirrer Couplings are not suitable for viscous media.

7. Waste disposal

The Globe Stirrer Coupling as well as its packing are made of recyclable materials. Please take further information on the materials used from the following specifications.

BOLA Globe Stirrer Couplings

8. Specifications

Bola Globe Stirrer Couplings

Cat.-No.	Opening for stirrer shaft mm	Upper dia. mm	Total length mm	Permissible speed rpm	Misalignment of axes mm	Transmittable torque Ncm	Material
C 394-01	Ø 4,0	Ø 8	190	1200	10	300	POM, glass-fibre reinforced PTFE, Stainless Steel, NBR (Perbunan)
C 394-02	Ø 6,5 und Ø 10,0	Ø 8	190				
C 394-03	Ø 8,0 und Ø 10,0	Ø 8	190				
C 398-04	Ø 4,0	Ø 10	190				
C 398-08	Ø 6,5 und Ø 10,0	Ø 10	190				
C 398-12	Ø 8,0 und Ø 10,0	Ø 10	190				
C 399-12	square 6	SW 8	180				
C 399-13	GL 10	Ø 10	170				

9. Spare parts / Accessories

» see page 25

Further accessories for optimum stirring and mixing can be found on our homepage www.bola.de.

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BOHLENDER GmbH