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Operating Instructions

Mode d'emploi

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BOLA Cordless Pumps for Acids and Caustig Solutions

BOLA Cordless Pumps for Acids and Caustic Solutions, Cat.-No.: G 870-01 and G 870-11.

OPERATING INSTRUCTIONS

| | |
|----------------------------------|----------|
| 1. Supplied with | 6 |
| 2. Security Advice | 6 |
| 3. Taking into operation | 6 |
| 4. Mode of Function | 7 |
| 5. Service / Maintenance | 7 |
| 6. Disturbances / Repairs | 7 |
| 7. Waste Disposal | 8 |
| 8. Specification | 8 |
| 9. Registered Trademarks | 8 |

Please read the information listed below thoroughly. We urgently ask you to observe the information about security, use and maintenance.

BOLA Cordless Pumps for Acids and Caustic Solutions can be used either with 2 batteries or accumulators Mono 1.5 V (batteries are not included in delivery). Due to a better capacity and for environmental reasons we recommend to use rechargeable batteries (= accumulators). BOLA Cordless Pumps for Acids and Caustic Solutions are available either with a suction pipe length of 400 mm or 600 mm.

BOLA Cordless Pumps for Acids and Caustic Solutions

1. Supplied with


- 1 BOLA Cordless Pumps for Acids and Caustic Solutions, optionally with suction pipe length 400 or 600 mm
- 1 Outlet Pipe
- 1 Operating Instructions

2. Security Advice

- » Do not take a damaged pump into operation.
- » The pump is suitable only for watery products without any solids.
- » Do not use the pump in explosive areas.
- » Do not pump products which charge electrostatically.
- » Do not use the pump without mounted outlet pipe.
- » After pumping, residues inside the pump can run out even if the pump is switched off.
- » The "OFF-switch" does not have any function. Upon letting the "ON-switch" off, the pump switches off.

3. Taking into operation


- » Remove the red battery housing (1) from the pump housing (2) by turning it counterclockwise.

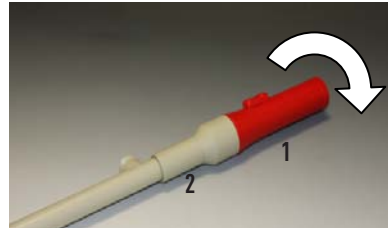
- » Fit two loaded batteries type Mono 1.5 V in the battery housing (1). The positive poles have to be in direction of the opening of the battery housing.

- » Screw the battery housing (1) on the pump housing (2) by turning it clockwise.

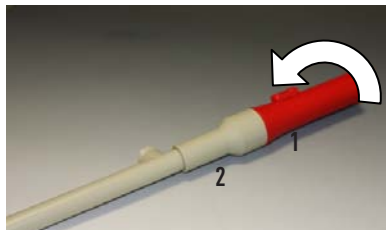
- » Mount the outlet pipe (3) by screwing the connecting nut (4) to the suction pipe (5).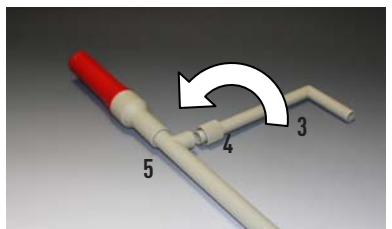

BOLA Cordless Pumps for Acids and Caustic Solutions

- » Introduce the suction pipe (5) into the liquid that has to be pumped. The end outlet pipe (3) has to be put into a collecting vessel (6).
- » Start to pump by pressing the "ON-button" (7). Once the "ON-button" is released, pumping is stopped immediately.

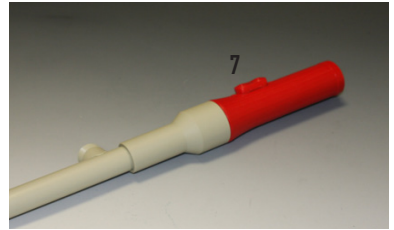4. Mode of Function

The pump is actuated by two standard batteries (accumulators) type Mono 1.5 V. The rotation of the pump shaft is transmitted by an encapsulated magnetic coupling to the pump impeller. This magnetic coupling protects at the same time the electronic components from being damaged by the pumped products.

5. Service / Maintenance

- » Rinse the pump with clear water after use.
- » Take off the accumulators if the pump is not used for a longer period of time.
- » Store the pump in suspended state.

6. Disturbances / Repairs

The following information shall help you to recognise and, if possible, to eliminate possible reasons for disturbances by yourself. If you are not able to handle the disturbance, please give us a call. We will then try to help you this way. If then, a solution to your problem cannot be found, we will ask you to return the cleaned pump in the original packing to us. Then we will check whether a cost-effective repair is possible.

BOLA Cordless Pumps for Acids and Caustig Solutions

| Objection | Reason and help |
|----------------------------------|--|
| The pump does not start to work. | <ul style="list-style-type: none"> » The batteries/accumulators are empty. Check whether the accumulators have to be charged and exchange them against new ones if necessary. » The batteries/accumulators are not inserted correctly. Check the direction of the poles. The positive poles have to be in direction of the opening of the battery housing (1). » The pump shaft is stuck. Rinse the pump with hot water. » The pump has sucked in solid particles. The pump is not suitable for liquids which contain solids. Rinse the pump with hot water. |
| The pump does not suck. | <ul style="list-style-type: none"> » The charge state of the batteries is too low. Check your accumulators and charge or exchange them if necessary. » The suction pipe is not deep enough immersed in the liquid. Immerse the suction pipe for at least 15 mm. » The product is too viscous. The pump is suitable only for watery products. |

7. Waste Disposal

BOLA Cordless Pumps for Acids and Caustic Solutions as well as their packing are made of recyclable materials. Please take further information on the materials used from the following specifications:

8. Specification

BOLA Cordless Pumps for Acids and Caustic Solutions

| Cat.-No. | Length of suction pipe | Suction pipe O.D. | Flow rate | Temperature resistance | Weight | Materials | Power Supply |
|----------|------------------------|-------------------|----------------------------------|------------------------|------------------|--------------------------------------|--|
| G 870-01 | 400 mm | approx. 25 mm | Max. 6 Liter pro Minute | + 5 °C bis + 60 °C | approx. 500 g | Polypropylene Hastelloy®, PTFE | Two accumulators, Type Mono 1,5 V |
| G 870-11 | 600 mm | | | | approx. 600 g | | |

9. Registered Trademarks

Hastelloy® is a registered trademark of Haynes International, Inc.

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BOHLENDER GmbH