Bedienungsanleitung
Operating Instructions
Mode d'emploi

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
OPERATING INSTRUCTIONS

1. Supplied with	7
2. Security instructions	7
3. Taking into operation	7
4. Mode of function	8
5. Service / Maintenance	8
6. Waste disposal	8
7. Customizations	8
8. Specifications	9

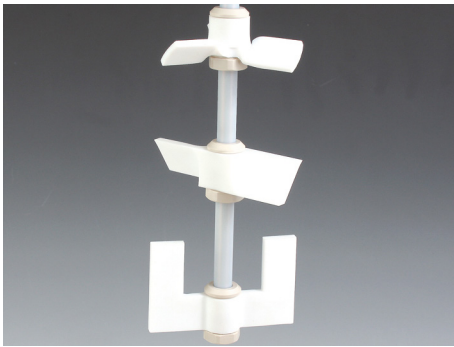Please read the information listed below thoroughly. We urgently ask you to observe the information about security, use and maintenance.

BOLA Blades

1. Supplied with

- 1 PTFE stirrer blade
- 1 Sleeve
- 1 Adjusting nut
- 1 Fork wrench
- 1 Operating instructions

2. Security instructions

- » Please do not take a damaged stirrer blade into operation.
- » Risk of injury from rotating pieces: all blades must be inside the stirring vessels during agitation.
- » The agitator must not be started before all rotating blades are inside the stirring vessel.
- » Only wear tight clothes and tight gloves to avoid being drawn in by rotating components. If you have long hair, please don't wear it loose.

3. Taking into operation

Please remove the adjusting nut (1) before assembling the blade on a stirrer shaft.




- 1 adjusting nut
- 2 sleeve
- 3 PTFE blade

Push the adjusting nut (1) on the stirrer shaft (4); hexagon downwards.

Then push the stirrer blade (3) with the sleeve (2) from above on the stirrer shaft (4), hexagon upwards. The hexagon of the sleeve (2) and the hexagon of the blade (3) can now be locked. Now screw the nut (1) and the sleeve (2) together with two rotations.


Note: If necessary for your application, you can assemble the complete stirrer blade overhead. The stirring direction however does not change.

Adjust the pre-assembled blade with the sleeve to the requested height.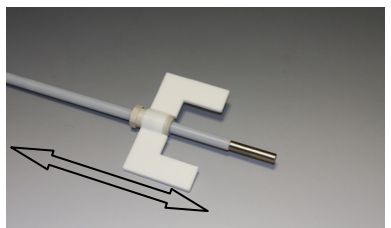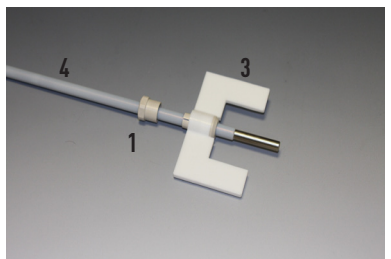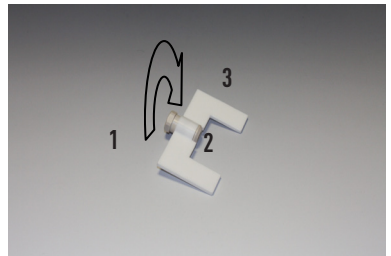


BOLA Blades


Screw up the adjusting nut (1) with the sleeve (2) in clockwise direction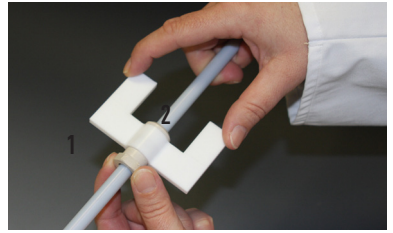

Tighten the adjusting nut (1) with the fork wrench (5). The adjusting nut is tight enough if there is no visible gap between nut and blade.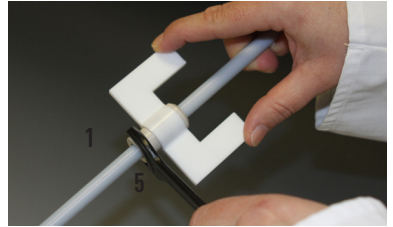

Attention: If you are using hard-walled stirrer shafts, for example made of glass or stainless steel, you must not flush the adjusting nut (1) with the blade. The nut must be fixed very carefully, and a gap of approx. 3 mm will remain visible between adjusting nut and blade. This gap does not have any influence on the stirring function.

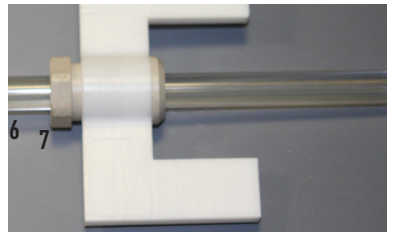4. Mode of function

The solid PTFE stirrer blades can be fixed on stirrer shafts with the indicated diameters. Twelve holding points provide the necessary clamping power.

5. Service / Maintenance

We recommend to detach the blade before cleaning. Please use commercial, non-abrasive detergents.

6. Waste disposal

The blades as well as the packing are made of recyclable materials. Please take further information about materials used from the following specifications.

7. Customizations

If you need a different version of a BOLA Stirrer Blade, do not hesitate to send us your inquiry including all requested features. As manufacturers, we are able to offer an individual custom manufacture.

In general, following modifications are possible:

- » a different stirring diameter
- » a different shaft diameter
- » a different material
- » a different stirring direction
- » a different shape of blade

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BOHLENDER GmbH